



Swift Normalase™ Amplicon Panels (SNAP) ***SARS-CoV-2, SARS-CoV-2 Additional Genome Coverage, and*** ***SARS-CoV-2 S Gene Panel***

Whole viral genome targeted next generation sequencing (NGS) assay

Single-tube multiplex PCR NGS library prep with high-throughput workflow

Protocol for Cat. Nos.:

COVG1-96, SARS-CoV-2 Panel (primers only, 96 rxns)

COVG1V2-96, SARS-CoV-2 Additional Genome Coverage Panel (primers only, 96 rxns)

SGENE-96, SARS-CoV-2 S Gene Panel (primers only, 96 rxns)

DNA Library Core Kits

Swift Normalase Amplicon Panel (SNAP) Core Kit (96 rxns)

Swift Normalase Amplicon Panel (SNAP) Core Kit (4x96 rxns bundle)

SNAP Indexing Primers

SNAP Set 1A to Set 2B Combinatorial Dual Indexing Primers (96-plex to 384-plex)

SNAP Unique Dual Indexing Primer Plates (96-plex to 1536-plex)

See Appendix Section H for details on Product Ordering Information.

Visit [swiftbiosci.com/protocols](https://www.swiftbiosci.com/protocols) for updates



Table of Contents

Product Information.....	2
Applications and Sample Types	2
Swift Normalase™ Amplicon Panels (SNAP) workflow.....	3
Kit Contents.....	4
Storage and Usage Recommendations.....	4
Materials and equipment not included.....	5
Reverse transcriptase and starting material recommendations.....	5
Tips and Techniques	5
Prepare SNAP libraries	6
Multiplex PCR Step.....	6
Panel-Specific Multiplex PCR Reaction Mix.....	6
Indexing PCR Step.....	7
Size selection and clean-up step 1	7
Size selection and clean-up step 2.....	8
Library Quantification	9
Introduction to Swift Normalase methods.....	9
Normalase I: Enzymatic Selection.....	10
Equal volume library pooling.....	10
Normalase II: Enzymatic Normalization	11
Normalase inactivation.....	12
Quantification and calibration of Swift Normalase pools.....	12
Appendix.....	13
Section A: Sequencing Recommendations.....	13
Section B: sequencing data analysis.....	13
Section C: Low Viral Load Input Recommendations.....	14
Section D: Indexing Primer Sequences	15
Section E: Swift 1536 UDIs - Plate Specification and Dimensions	15
Section F: Plate Usage Guidelines	15
Section G: Plates Layout	16
Section H: Product Ordering Information.....	19
Revision History.....	20
Notice to Purchaser: Limited License	20

Support

For additional support please contact Swift at TechSupport@swiftbio.com, or by phone: 734.330.2568 (9:00 am – 5:00 pm ET, Monday through Friday).

Product Information

The SNAP SARS-CoV-2 Kits offer a robust NGS workflow that provides optimal coverage and NGS data quality on Illumina® sequencing platforms. These kits leverage patented multiplex PCR technology, enabling library construction from first- or second-strand cDNA using tiled primer pairs to target the 29.9 kb viral genome or S gene along with a single pool of multiplexed primer pairs. Primers were designed against the NCBI Reference Sequence NC_045512.2 (Severe acute respiratory syndrome coronavirus 2 isolate Wuhan-Hu-1, complete genome).

Features	Specifications
Design coverage and panel information	SARS-CoV-2 Panel 98.0% (29,313 of 29,903 total bases) 341 amplicons, sized 116-255 bp (average 150 bp)
	SARS-CoV-2 Additional Genome Coverage Panel 99.7% (29,828 of 29,903 total bases) 345 amplicons, sized 116-255 bp (average 150 bp)
	SARS-CoV-2 S gene Panel 100% of S gene (100% of S gene bases) 43 amplicons, sized 140-150 bp (average 146 bp)
Input Material	1 st or 2 nd strand cDNA Minimum 10 – 100+ viral copies (RT-qPCR Ct value 30-40)
Time	2 hours cDNA-to-Library
	3 hours cDNA-to-Normalized-Library Pool
Components	Included: <ul style="list-style-type: none"> • Target-specific multiplex primer pool • PCR and library prep reagents Optional: <ul style="list-style-type: none"> • Combinatorial Dual Indexed Adapters • Swift Normalase Note: Kits do not include RT module or magnetic beads
Multiplexing Capability	Up to 384 CDIs or 1536 UDIs
Recommended Depth	250-500k reads/library for SARS-CoV-2 and SARS-CoV-2 Additional Genome Coverage Panels 0.25M reads/library for SARS-CoV-2 S Gene Panel

*Please inquire with your Swift sales representative or distributor to obtain a copy of the primer design file.

Applications and Sample Types


- **Applications:** detection, variant calling, epidemiological studies, public health surveillance, among others.
- **Internally Demonstrated Sample Types:** nasopharyngeal/oropharyngeal swabs, sputum, bronchoalveolar lavage (BAL), stool, and wastewater, among others.

Swift Normalase™ Amplicon Panels (SNAP) workflow

Swift Normalase Amplicon Panels (SNAP) utilize multiple overlapping amplicons in a single tube, using a rapid, 2-hour workflow to prepare ready-to-sequence libraries. The PCR1+PCR2 workflow generates robust libraries, even from 10–100 viral genome copies. The libraries may be quantified with conventional methods such as Qubit® or Agilent Bioanalyzer and normalized by manual pooling or normalized enzymatically with the included Swift Normalase reagents.

This protocol includes instructions for a Multiplex PCR step to enrich target sequences, an Indexing PCR step to amplify and add combinatorial or unique dual indexed adapter sequences, and an optional downstream Swift Normalase step to produce an equimolar library pool.

SNAP Workflow

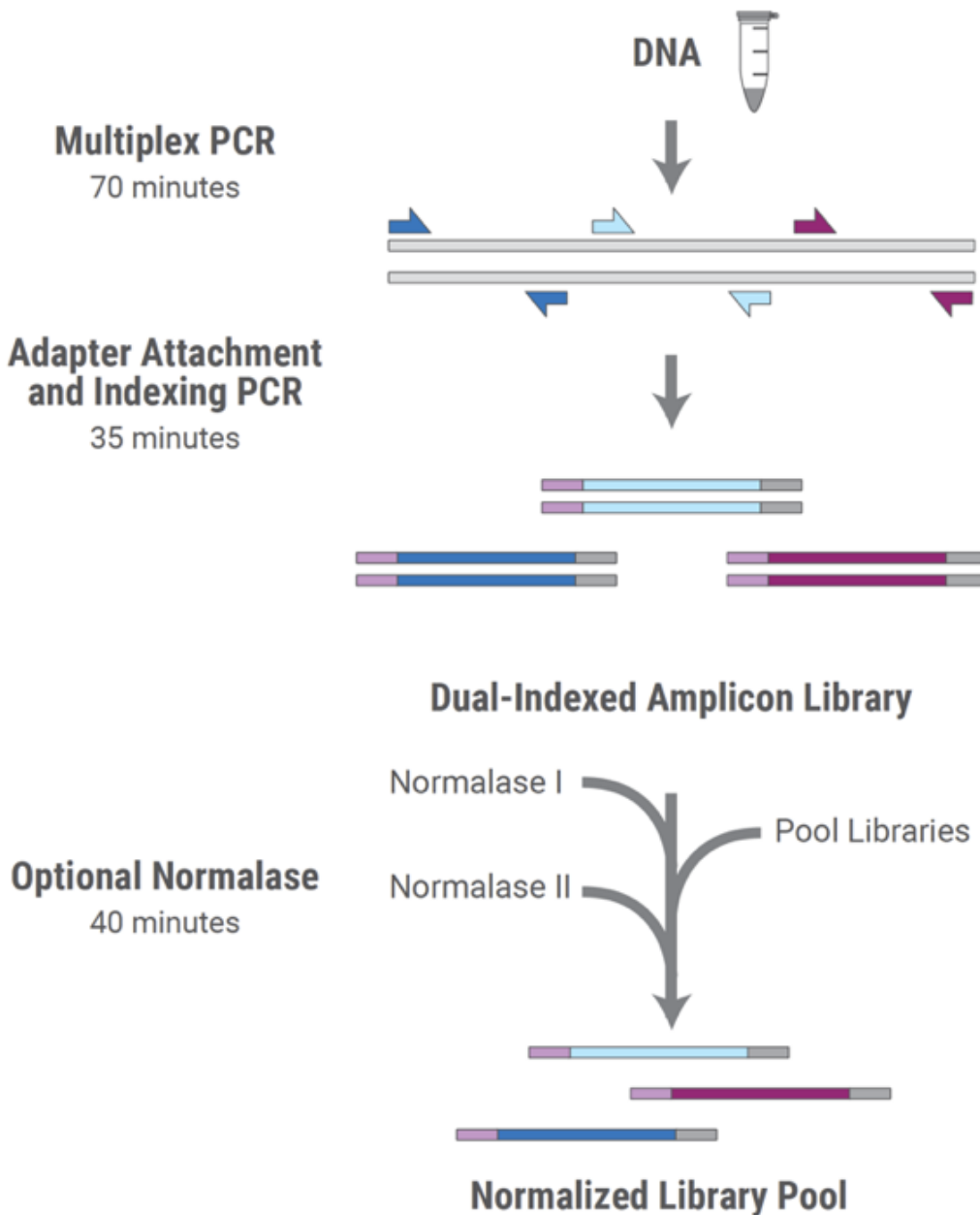

Kit Contents

The SNAP Kits contain sufficient reagents for the preparation of 96 libraries (10% excess volume provided).

Protocol Stage	Component	96 rxns	Storage
Multiplex PCR (Pre-PCR box)	● Reagent G1*	211 µL	-20°C***
	● Reagent G2	317 µL	
	● Enzyme G3	1,584 µL	
	● Pre-PCR TE	1,200 µL	
Indexing PCR (Post-PCR box)	● Reagent I1	348 µL	
	● Enzyme I2	53 µL	
	● Enzyme I3	15 µL	
	● Enzyme I4	2,640 µL	
	● Post-PCR TE	4,161 µL	
Indexing Primer Box**	● SNAP CDI D50X	26 µL each	
	● SNAP CDI S7XX	15 µL each	
Indexing Primer Plates**	● SNAP UDI (pre-mixed pairs)	12 µL per well	
Normalase (Post-PCR box)	● Buffer S1	454 µL	
	● Reagent S2	21 µL	
	● Enzyme S3	53 µL	
	● Buffer N1	506 µL	
	● Enzyme N2	21 µL	
	● Reagent X1	105 µL	
Additional Reagents	PEG NaCl	20 mL	Room Temp

*Reagent G1 is included in Cat. No. COVG1-96, COVG1V2-96, and SGENE-96. Additional Pre-PCR and Post-PCR box reagents and PEG NaCl are included in Cat. Nos. SN-5X296 and SN-5X384.

** Indexing boxes are Cat. Nos. SN-5S1A96, SN-5S1B96, SN-5S2A96 and SN-5S2B96. Indexing Primer Plates are Cat. Nos. SN91096-1-PLATE, SN91096-2-PLATE, SN91096-3-PLATE, SN91096-4-PLATE (available as individual plates), and SN91384-PLATES, SN91384-B-PLATES, SN91384-C-PLATES and SN91384-D-PLATES (available as 4x96 bundle plates).

***Temperature range -15°C to -25°C.

Storage and Usage Recommendations

- Upon receipt, store the SNAP Library Kit products at -20°C with the exception of the PEG solution, which is stored at room temperature.
- Separate the multiplex PCR Reagents (keep in pre-PCR area) from the Indexing and Swift Normalase Reagents (keep in post-PCR area).
- To maximize use of enzyme reagents when ready to use, remove enzyme tubes from -20°C storage and place on ice for 10 minutes prior to pipetting. Attempting to pipette enzymes at -20°C may result in shortage of enzyme reagents.
- After thawing reagents on ice, briefly vortex (except enzymes) to mix well, then pulse spin to collect contents before proceeding. Enzyme G3 is the only enzyme that may be vortexed.
- Always add reagents to the master mix in the specified order as stated throughout the Protocol. The indexing primers (SNAP CD and UD Indexes) are the only reagents that are added individually to each sample.

Materials and equipment not included

- Reverse Transcriptase (RT) Module—See below
- SPRIselect beads (Beckman Coulter, Cat. Nos. B23317/B23318/B23319) or Agencourt AMPure XP beads (Beckman Coulter, Cat. Nos. A63880/A63881/A63882)
- Permagen Magnetic Separator (Cat. No. MSR812), Agencourt SPRIPlate were used for internal benchmarking. Similar magnetic rack for magnetic bead clean-ups can be substituted.
- Qubit®, Nanodrop, or other similar input RNA quantification assay
- qPCR-, electrophoretic-, or fluorometric-based library quantification assay for Illumina® libraries
- Microcentrifuge
- Vortex
- Programmable thermocycler
- 0.2 mL PCR tubes or 96-well plates
- Aerosol-resistant tips and pipettes ranging from 1 to 1000 µL
- 200 proof (absolute) ethanol and nuclease-free water (both molecular biology grade) for preparation of fresh 80% ethanol
- Pipette tips (e.g., 8-channel or 12-channel), 8-tube strips, an un-skirted 96 well plate, or plate puncher for pre-piercing the foil seal if using single-use UD indexing plates.

Reverse transcriptase and starting material recommendations

The Swift SARS-CoV-2 amplicon panels support first- or second-strand cDNA as input. Select a cDNA synthesis module that supports two-step RT-PCR, includes random primers, and has a processivity of >1 kb.

The amplicon panels have been demonstrated with synthetic SARS-CoV-2 genome (Twist Biosciences Cat. No. 102024). As few as 10–100 viral genome copies were mixed with 50 ng of Universal Human Reference RNA (Agilent 740000) and used as input into the Superscript® IV First-Strand Synthesis System (Thermo Fisher Cat. No. 18091050). The manufacturer's protocol was followed as written using the random primers and associated specifications, with a single modification of increasing the RT incubation step (50°C) from 10 min to a minimum of 30 min. Optional host gDNA/RNA removal steps were not performed. The optional RNase H step was not performed, and 10 µL of the resulting cDNA was used directly as input into the amplicon panel.

Tips and Techniques

Assemble all reagent master mixes and reactions ON ICE and scale volumes as appropriate, using 5% excess volume to compensate for pipetting loss. Calculate the total volume of the master mixes and prepare them in advance to ensure the magnetic beads do not over-dry during size selection steps while awaiting completion of master mix assembly. Neglecting to store master mixes and reagents on ice prior to incubations reduces yields and performance of this product.

This workflow, like any amplicon enrichment technology, poses a risk of contamination of surfaces and other samples following the amplification step. Please use caution when opening your sample tubes following the Multiplex PCR step. To reduce the risk of RNA and library contamination, physically separate the laboratory space, equipment, and supplies where pre-PCR and post-PCR processes are performed. Follow the instructions below to avoid cross-contamination:

- Clean lab areas using 0.5% sodium hypochlorite (10% bleach).
- Use barrier pipette tips to avoid exposure to potential contaminants.
- Always change tips between each sample.

Prepare SNAP libraries

Multiplex PCR Step

1. Load the Multiplex PCR program and allow the block to reach 98°C before loading samples (**confirm lid heating is turned ON and is set to reach 105°C**).

Pre-Program Thermocycler		
	98°C, 30 sec	
Multiplex PCR	98°C, 10 sec	4 cycles
	61°C, 5 min	
	65°C, 1 min	
	98°C, 10 sec	
	64°C, 1 min	18 cycles*
	65°C, 1 min	
	4°C, Hold	

* The PCR cycle number can be increased for samples that may give low yields. See the [Low viral load input recommendations](#) for adjusted thermocycling programs (Appendix, Section C).

2. Gently nutate Enzyme G3 at Room Temperature for 5 minutes or until all solutes appear to be in solution. Place back on ice for remainder of use.
3. Load 10 µL of cDNA sample directly into each PCR tube.
4. Keep all tubes on ice during assembly of the master mix and the reaction until placed in thermocycler.

Panel-Specific Multiplex PCR Reaction Mix

Before mixing, calculate the total volume of the master mix based on the number of reactions required with appropriate overage for pipetting. Vortex components G1 and G2 and pulse-spin tubes to collect contents. Make the Multiplex PCR Reaction Mix as shown in the below table. Keep prepared master mix on ice until ready to use.

Reagents	Volume per Sample
● Reagent G1*	2 µL
● Reagent G2	3 µL
● Enzyme G3	15 µL
Reaction Mix	20 µL

* Reagent G1 is the panel-specific set of amplification primers.

5. Mix the master mix well and then add 20 µL of the Multiplex PCR Reaction Mix to each 10 µL input cDNA sample on ice. Mix well, then place in the thermocycler and run the program.

IMPORTANT: Move samples to post-PCR area before opening tubes.

6. Near the completion of the thermocycler run, prepare the Indexing Reaction Mix (see below) in the post-PCR area. Assemble this reaction mix on ice and keep it cold until adding it to samples in the Indexing Step. All components except indexes should be master-mixed when running multiple samples in parallel.

Indexing PCR Step

Before mixing, calculate the total volume of the master mix based on the number of reactions, with appropriate overage for pipetting. We recommend preparing at least 10 reactions at any one time to maintain a volume of Enzyme I3 that can be accurately pipetted. Keep prepared master mix on ice.

Reagents	Volume per Sample
• Reagent I1	3.3 μ L
• Enzyme I2	0.5 μ L
• Enzyme I3	0.1 μ L
• Enzyme I4	25 μ L
Reaction Mix	28.9 μL

IMPORTANT: Keep prepared master mix ON ICE during size selection and clean-up step 1.

Size selection and clean-up step 1

7. Ensure beads and samples are at room temperature. Briefly vortex beads to homogenize before use.
8. Add **30 μ L (ratio: 1.0)** of magnetic beads to each 30 μ L sample. Mix by vortexing. Pulse-spin the samples in a microfuge. Ensure no bead-sample suspension droplets are left on the sides of the tube.
9. Incubate the samples for 5 minutes at room temperature off the magnet.
10. Place the sample tubes on a magnetic rack until the solution clears and a pellet is formed (~5 minutes).
11. While leaving your sample on the magnet, remove and discard the supernatant without disturbing the pellet (approximately 5 μ L may be left behind). Leave tubes on the magnet.
12. Add 180 μ L of freshly prepared 80% ethanol solution to the pellet while it is still on the magnet. Use care not to disturb the pellet. Incubate for 30 seconds, and then carefully remove the ethanol solution.
13. Repeat, for a second wash with the ethanol solution.
14. Pulse-spin the samples in a microfuge, place back onto the magnet, and remove any residual ethanol solution from the bottom of the tube with a small volume tip.
15. Resuspend each bead pellet in 17.4 μ L Post-PCR TE Buffer. Proceed to the Indexing PCR step.

IMPORTANT: Continue working in the post-PCR area. Keep samples at room temperature. At no time should “with bead” samples be stored on ice, as this affects binding to magnetic beads.

16. Load the Indexing PCR program and allow the block to reach 37°C before loading samples (**confirm lid heating is turned ON and is set to reach 105°C**).

Thermocycler Program		
	37°C, 20 min	
	98°C, 30 sec	
Indexing PCR	98°C, 10 sec	9 cycles*
	60°C, 30 sec	
	66°C, 1 min	
	4°C, Hold	

* The PCR cycle number can be increased for samples that may give low yields. See the [Low viral load input recommendations](#) for adjusted thermocycling programs (See Appendix, Section C).

17. Add a unique combination of 2 µL SNAP CD Index D50X + 1.7 µL SNAP CD Index S7XX to each sample, OR
Add 3.7 µL of a pre-mixed SNAP UD Indexing primer pair to each sample if using the single use plates (See [Appendix, Section F for UDI plate usage guidelines](#)).
18. Add 28.9 µL of the cold Indexing PCR Reaction Mix to each sample and mix thoroughly (total volume 50 µL).
19. Place in the thermocycler and run the program.

Size selection and clean-up step 2

20. Ensure PEG NaCl solution is at room temperature. Briefly vortex the PEG NaCl solution to homogenize before use.
21. Add 32.5 µL (ratio: 0.65) of PEG NaCl solution to each 50 µL sample. Mix by vortexing. Ensure no bead-sample suspension droplets are left on the sides of the tube.
22. Incubate the samples for 5 minutes at room temperature off the magnet.
23. Pulse-spin the samples in a microfuge. Place the sample tubes on a magnetic rack until the solution clears and a pellet is formed (~5 minutes).
24. While leaving your sample on the magnet, remove and discard the supernatant without disturbing the pellet (approximately 5 µL may be left behind). Leave tubes on the magnet.
25. Add 180 µL of freshly prepared ethanol solution to the pellet while it is still on the magnet. Use care not to disturb the pellet. Incubate for 30 seconds and then carefully remove the ethanol solution.
26. Repeat step 25 for a second wash with the ethanol solution.
27. Pulse-spin the samples in a microfuge, place back onto the magnet, and remove any residual ethanol solution from the bottom of the tube with a small volume tip.
28. Proceed immediately to add 20 µL of Post-PCR TE buffer and resuspend the pellet, mixing well by pipetting up and down until homogenous. Incubate at room temperature for 2 minutes off the magnet. Then place the sample back on the magnet and transfer the clean 20 µL library eluate to a fresh tube. Ensure that eluate

does not contain magnetic beads (indicated by brown coloration in eluate). If magnetic beads are present, pipette eluate into a new tube, place on magnet, and transfer eluate again.

NOTE: Libraries are now completed and ready to sequence. Please proceed with quantifying and pooling libraries, using either conventional fluorometric (e.g., Qubit) or electrophoretic (e.g., Bioanalyzer) methods, or proceed to Normalase below (reagents are included).

Library Quantification

Accurate library quantification is essential to load the sequencing instrument properly. Libraries can be quantified using fluorometric, electrophoretic, or qPCR-based methods and normalized manually. Alternatively, libraries can be enzymatically normalized following the Swift Normalase protocol below. Note, for optimal normalization using the Swift Normalase method, a minimum of 12 nM yield is needed per sample. If library yields are below 12 nM, increase the number of PCR cycles to pass the 12 nM threshold or switch to the 6 nM threshold Swift Normalase protocol described below.


Introduction to Swift Normalase methods

Here we provide instructions for optional enzymatic normalization of multiplexed SNAP NGS libraries for equimolar pools and balanced sample representation in sequencing. The protocol is designed for SNAP libraries that produce consistent amplified library yields of ≥ 12 nM following Indexing PCR, and it generates an equimolar library pool. Most samples processed with SNAP protocol produce amplified library yields of 12 nM or greater; however, if there is concern that not all samples will reach 12 nM, adjusting Swift Normalase chemistry to require a minimum of only 6 nM can alternatively be performed.

A simple [calculator](#) for converting between ng/ μ L and nM is available. For the SNAP SARS-CoV-2 panels, please use a finished library size of 285 bp (for use in the Base Pair Length column).

The workflow consists of three steps for libraries amplified to a minimum yield of 12 nM during the adapter attachment and Indexing PCR step:

1. **Normalase I** to select a 4 nM (or 2 nM if using ≥ 6 nM option) library fraction enzymatically
2. **Library Pooling** of samples for multiplexed sequencing
3. **Normalase II** to generate an equimolar library pool enzymatically

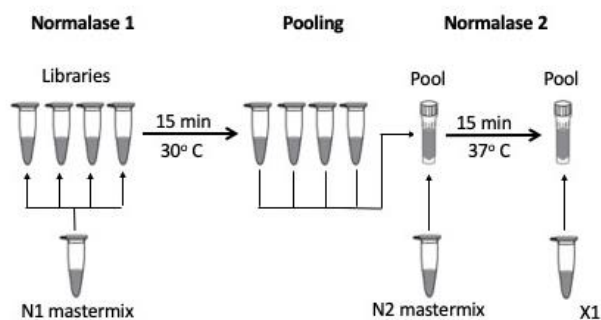

Workflow schematic: Normalase I master mix is added to samples and incubated at 30°C for 15 minutes. Sample pooling is then performed, and then the Normalase II master mix is added to the pool. Samples are incubated at 37°C for 15 minutes. Reagent X1 inactivates the reaction, and a final equimolar pool is produced.

Notes regarding Swift Normalase specification

The Swift Normalase product specification is defined by cluster density of the Swift Normalase pool when loaded on a MiSeq v2 flow cell at 12 nM to achieve a 1,000–1,200 K/mm² cluster density and CV $\leq 15\%$ within a pool. Across Illumina platforms, library types, and insert sizes, the optimization of loading concentration may be required to achieve the optimal number of reads supported by the flow cell of choice.

Normalase I: Enzymatic Selection

If you are concerned that the 12 nM threshold has not been met after Indexing PCR:

- Spot check library yields using either fluorometric methods (e.g., Qubit) or electrophoretic methods (e.g., Bioanalyzer).
 - A Swift Normalase workflow modification can be performed that requires a 6 nM threshold to obtain a 2 nM Swift Normalase pool (see below).
1. Pre-set a thermocycler program as listed below.

Thermocycler Program
15 min at 30°C with open lid or lid heating OFF

2. Prepare the Normalase I master mix as listed in the table below. The mix can be prepared at room temperature and stored on ice until use if prepared in advance. Ensure that it is thoroughly mixed by moderate vortexing followed by a pulse spin to collect contents prior to use. For libraries with lower yields \geq 6 nM, or for a final pool of 2 nM (instead of 4 nM), please use half of the specified volume of Reagent S2 and add an equal volume of TE, thus reducing concentration by two-fold, then proceed as written.

Reagent	Per library	24 Libraries	96 Libraries
• Buffer S1	4.3 μ L	103.2 μ L	412.8 μ L
• Reagent S2	0.2 μ L	4.8 μ L	19.2 μ L
• Enzyme S3	0.5 μ L	12 μ L	48 μ L
Total volume	5 μL	120 μL	480 μL

IMPORTANT: The Normalase I master mix should be built for a minimum of 10 reactions to ensure pipetting accuracy.

3. Using a calibrated P10 pipette, add 5 μ L of Normalase I master mix into each 20 μ L library eluate at room temperature and thoroughly mix by moderate vortexing for 5 seconds.
4. Spin down the sample tube in a microfuge. Place in the thermocycler and run the program.

----- **SAFE STOPPING POINT** -----

Store freshly prepared libraries at -20°C post-Normalase I.

Equal volume library pooling

Sufficient Normalase II reagents are supplied so this step can be repeated to enable various re-pooling combinations, as only 5 μ L of post-Normalase I library (out of a 25 μ L volume) is used for pooling. Also note that stability of normalized pools (after Normalase II) is limited with a storage time of four weeks, since the resulting normalized pools contain single-stranded DNA. Therefore, if re-sequencing is required after four weeks, for best results re-pool the Normalase I libraries and repeat Normalase II and inactivation.

NOTE: If you are pooling fewer than five libraries, contact tech support at TechSupport@swiftbio.com for low-plex pooling recommendations.

NOTE: If pooling 5 μL per sample does not generate a normalized pool of sufficient volume for instrument loading, contact tech support at TechSupport@swiftbio.com for high sample volume pooling recommendations.

IMPORTANT: Consider your desired number of reads for each sample and pool only those samples together that have the same required depth. For example, samples each requiring 50,000 reads can be pooled together, whereas samples requiring 1 million reads should be combined in a separate pool. Thus, you can adjust your ratio of pools when loading the instrument to achieve the desired sequence depth for each pool.

1. Following the Normalase I incubation, generate a library pool (or pools) by placing 5 μL of each individual library into a 0.2 mL PCR tube if pooling 30 libraries or less (this achieves a final volume of up to 186 μL). Alternatively, use a 1.5 mL screw-cap microfuge tube, particularly when pooling more than 30 libraries, as the volume will exceed the PCR tube maximum volume.

To ensure even pooling, use of a calibrated P10 pipette will produce the best results.

2. Thoroughly mix, spin the library pools in a microfuge, and proceed to the Normalase II reaction.

Normalase II: Enzymatic Normalization

1. Pre-set a thermocycler program as listed below. Alternatively, if using a 1.5 mL screw cap microfuge tube, set a heat block at 37°C.

Thermocycler Program	Heat Block (1.5 mL screw cap microfuge tube)
15 min at 37°C with open lid or lid heating OFF	15 min at 37°C

IMPORTANT: Pre-mix Normalase II master mix (listed in the table below). The master mix can be stored on ice until use, and then added to pools at room temperature.

Reagents*	Per Library	24 Libraries	96 Libraries
• Buffer N1	0.96 μL	23.04 μL	92.16 μL
• Enzyme N2	0.04 μL	0.96 μL	3.84 μL
Total Volume	1 μL	24 μL	96 μL

*It is recommended to prepare Normalase II master mix for 24 samples even if you are processing less than 24 samples in order to avoid pipetting extremely low volumes; for best results use a calibrated P2 pipette for adding Enzyme N2. Although sufficient reagents are supplied for up to five repeated Normalase II reactions per sample, repeatedly processing a lower number of samples will result in significant loss of Normalase II reagents.

1. Add 1 μL of Normalase II master mix multiplied by the total number of libraries within each prepared pool.
2. Mix well by vortexing for 5 seconds and spin down the library pools in a microfuge.
3. Place the library pools in the thermocycler and run the program or place the 1.5 mL screw cap microfuge tubes into the 37°C heat block.

Normalase inactivation

1. Following the Normalase II reaction, pre-set a thermocycler program as listed below.

Thermocycler Program	Heat Block (1.5 mL screw cap microfuge tube)
Hold at 95°C 2 min at 95°C with lid kept at 95°C Hold at 4°C	2 min at 95°C

2. Add 0.2 µL of Reagent X1 multiplied by the total number of libraries within each prepared pool (see examples below).

Reagent	Per library	24-Plex pool	96-Plex pool
• Reagent X1	0.2 µL	4.8 µL	19.2 µL

3. Place the library pools in the thermocycler and advance the program or place the 1.5 mL screw cap microfuge tubes into the heat block. If using a 1.5 mL screw cap microfuge tube, set a heat block at 95°C to incubate your library pools, being careful not to incubate the samples longer than 2 minutes.
4. Your final multiplexed library pools are now equimolar. Proceed to qPCR quantification of your Swift Normalase pool and sequencing. It is not necessary to perform an additional purification step.

Quantification and calibration of Swift Normalase pools

To ensure optimal sequencing results, perform a qPCR quantification on your final Swift Normalase pool(s). Final library pools are ssDNA and cannot be quantified by dsDNA-based fluorometric methods or fragment analysis. If you do not have a qPCR assay, confirm a commercially available kit by calibrating your qPCR results and sequencer loading concentrations before proceeding (for example KAPA Library Quantification Kit).

The Normalase 4 nM formulation may not conform to your qPCR assay quantification due to the lack of precision across different qPCR assays and laboratory practices. Using a qPCR assay that reproducibly predicts an optimal number of reads on your sequencing instrument, load your final pool based on your qPCR results. Across other Illumina platforms, library types, and insert sizes, optimization of loading concentration may be required to achieve the optimal number of reads supported by the flow cell of choice. If you have chosen the 6 nM to 2 nM option but require a higher pool concentration for your sequencer, perform a 2.0X SPRI to concentrate pools and then proceed to qPCR quantification and loading.

Appendix

Section A: Sequencing Recommendations

SNAP libraries may be sequenced using paired-end sequencing on Illumina instruments. We recommend using 2 x 150 paired-end reads. The depth of coverage required will depend on the application. Be sure to use either 8 (CDI) or 10 (UDI) sequencing cycles for each index read.

For the SARS-CoV-2 Panel, >50,000 reads per library are recommended for presence/absence (+/-) confirmation, whereas 250–500k reads per library are recommended for variant calling. Accordingly, the following table shows examples of the number of libraries that can be multiplexed to achieve this depth per sequencing run:

		Libraries Per Sequencing Run					
		MiSeq®				MiniSeq®	
Panel	Application	v2 Nano	v2 Micro	v2	v3	Mid-Output	High-Output
SARS-CoV-2 Panels	+/- Confirmation	40	160	600	1,000	320	1,000
	Variant Calling	8	32	120	200	64	120
S gene	Variant Calling	40	160	600	1,000	320	600

Due to the complexity of the libraries, no PhiX spike-in is required on MiSeq or MiniSeq instruments. The NextSeq550 may be sensitive to low complexity, and PhiX or another suitable high-complexity library spike-in may be required. Contact Illumina technical support for further information regarding sequencing instrument compatibility with low-complexity sequences.

Section B: sequencing data analysis

Dockerized data analysis workflow

For customers using the SNAP SARS-CoV-2 Panel Kits who are comfortable with command line tools and want a ready-to-use variant calling analysis workflow which they can run on a local Linux machine, Swift offers a full variant calling workflow with all tools and reference files pre-installed and configured in a Docker image available for download.

Features

- One easy install: Docker image file can be downloaded and loaded onto a Linux server running the Docker daemon
- Easy-to-run analysis workflow includes a run script which handles all Docker commands, requiring only the input FASTQ files and the panel master file as arguments

Requirements

- Linux server with Docker installed
- At least 8 GB of RAM (≥ 32 GB recommended)

Analysis Overview

1. Adapter trimming and filtering out of reads < 30 bases long (Trimmomatic)
2. Alignment (BWA MEM)
3. Primer trimming (Primerclip)
4. On-target and coverage metrics calculation and reporting

5. Variant calling (GATK Haplotype Caller)

Follow this [Download Link](#) for the analysis workflow.

For more detailed information about primer trimming, please review the [Primerclip Technical Note](#).

Section C: Low Viral Load Input Recommendations

If your samples have a low viral load, please use the following adjustments to cycle number for both the Multiplex PCR and Indexing PCR thermocycling programs, as shown below in **blue underlined text**.

Multiplex PCR Step

1. Load the Multiplex PCR program and allow the block to reach 98°C before loading samples (**confirm lid heating is turned ON and is set to reach 105°C**).

Pre-program thermocycler		
	98°C, 30 sec	
Multiplex PCR	98°C, 10 sec	4 cycles
	61°C, 5 min	
	65°C, 1 min	
	98°C, 10 sec	
	64°C, 1 min	<u>24 cycles</u>
	65°C, 1 min	
	4°C, Hold	

Then proceed as written.

Indexing PCR Step

16. Load the Indexing PCR program and allow the block to reach 37°C before loading samples (**confirm lid heating is turned ON and is set to reach 105°C**).

Thermocycler program		
	37°C, 20 min	
Indexing PCR	98°C, 30 sec	
	98°C, 10 sec	<u>5 cycles</u>
	60°C, 30 sec	
	66°C, 1 min	
	4°C, Hold	

Then proceed as written.

Section D: Indexing Primer Sequences

The full-length indexing primer sequences are below, where the underlined text indicates the location of the index sequences, which are 8b for CDI and 10b for UDI. These sequences represent the adapter sequences following completion of the **Indexing PCR** step.

Index 1 (i7) Adapter:

5' – GATCGGAAGAGCACACGTCTGAACTCCAGTCACXXXXXXXXXX(XX)ATCTCGTATGCCGTCTTCTGCTTG – 3'

Index 2 (i5) Adapter:

5' – AATGATACGGCGACCACCGAGATCTACACYYYYYYYYY(Y)ACACTCTTCCCTACACGACGCTCTTCCGATCT – 3'

Please refer to the accompanying **Swift Index Master List** available at <https://swiftbiosci.com/protocols/> for index sequences for preparing your Illumina sequencing sample sheet on your instrument of choice. Contact TechSupport@swiftbio.com if you would like assistance confirming compatibility of your own primers with the SNAP workflow, or contact your local sales representative or distributor to inquire about the purchase of custom Swift Normalase Indexing Primers that use your own index sequences.

Section E: Swift 1536 UDIs - Plate Specification and Dimensions

This product is dispensed in a 96-well plate. Physical specifications are below:

<u>Plate Dimension</u>	<u>Low-Profile 96-Well Skirted Plates</u>
Length at base plane	127.76 mm
Width at base plane	85.48 mm
Height overall	16.06 mm
Well depth	14.81 mm
Well diameter at opening	5.46 mm
Well diameter at bottom of conical section	2.64 mm
Well volume	200 µL
Well spacing	9.00 mm
Well angle	17.5°
<i>Well offset</i>	
Left edge to well A1	14.38 mm
Top edge to well A1	11.24 mm
Left edge to H12	113.38 mm
Top edge to H12	74.24 mm

Section F: Plate Usage Guidelines

Prior to piercing the foil and pipetting out the necessary indexes, please be sure to thaw the plate at room temperature, vortex briefly, and centrifuge for one minute to spin down the primer reagents to the bottom of the plate wells.

Carefully pre-pierce the foil seal for the intended well(s) prior to pipetting the primer mix out of the plate to add to your reaction(s). Pre-piercing the foil avoids accidental clogging of pipette tips used for liquid pipetting and also avoids the introduction of foil into the reaction. In addition, pre-piercing the foil reduces the resistance to multi-channel pipettors which can result in undesired movement of the plate that may cause cross-contamination of

reagents. The foil may be pre-pierced with pipette tips (e.g., 8-channel or 12-channel), 8-tube strips, an un-skirted 96 well plate or plate puncher.

During the **Indexing PCR** step, use 3.7 μ L of a unique indexing primer pair (SU001-SU1536 UDIs) to amplify and index each library, where the UDI primer pair must be added individually to each sample.

Libraries made with uniquely indexed adapters may be pooled prior to cluster generation, subjected to the Swift Normalase chemistry, and co-sequenced on the same Illumina flow cell.

Section G: Plates Layout

SN91384-PLATES (All four plates; SN91096-1-PLATE, SN91096-2-PLATE, SN91096-3-PLATE, SN91096-4-PLATE are included in bundle).

SN91096-1-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU001	SU009	SU017	SU025	SU033	SU041	SU049	SU057	SU065	SU073	SU081	SU089
B	SU002	SU010	SU018	SU026	SU034	SU042	SU050	SU058	SU066	SU074	SU082	SU090
C	SU003	SU011	SU019	SU027	SU035	SU043	SU051	SU059	SU067	SU075	SU083	SU091
D	SU004	SU012	SU020	SU028	SU036	SU044	SU052	SU060	SU068	SU076	SU084	SU092
E	SU005	SU013	SU021	SU029	SU037	SU045	SU053	SU061	SU069	SU077	SU085	SU093
F	SU006	SU014	SU022	SU030	SU038	SU046	SU054	SU062	SU070	SU078	SU086	SU094
G	SU007	SU015	SU023	SU031	SU039	SU047	SU055	SU063	SU071	SU079	SU087	SU095
H	SU008	SU016	SU024	SU032	SU040	SU048	SU056	SU064	SU072	SU080	SU088	SU096

SN91096-2-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU097	SU105	SU113	SU121	SU129	SU137	SU145	SU153	SU161	SU169	SU177	SU185
B	SU098	SU106	SU114	SU122	SU130	SU138	SU146	SU154	SU162	SU170	SU178	SU186
C	SU099	SU107	SU115	SU123	SU131	SU139	SU147	SU155	SU163	SU171	SU179	SU187
D	SU100	SU108	SU116	SU124	SU132	SU140	SU148	SU156	SU164	SU172	SU180	SU188
E	SU101	SU109	SU117	SU125	SU133	SU141	SU149	SU157	SU165	SU173	SU181	SU189
F	SU102	SU110	SU118	SU126	SU134	SU142	SU150	SU158	SU166	SU174	SU182	SU190
G	SU103	SU111	SU119	SU127	SU135	SU143	SU151	SU159	SU167	SU175	SU183	SU191
H	SU104	SU112	SU120	SU128	SU136	SU144	SU152	SU160	SU168	SU176	SU184	SU192

SN91096-3-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU193	SU201	SU209	SU217	SU225	SU233	SU241	SU249	SU257	SU265	SU273	SU281
B	SU194	SU202	SU210	SU218	SU226	SU234	SU242	SU250	SU258	SU266	SU274	SU282
C	SU195	SU203	SU211	SU219	SU227	SU235	SU243	SU251	SU259	SU267	SU275	SU283
D	SU196	SU204	SU212	SU220	SU228	SU236	SU244	SU252	SU260	SU268	SU276	SU284
E	SU197	SU205	SU213	SU221	SU229	SU237	SU245	SU253	SU261	SU269	SU277	SU285
F	SU198	SU206	SU214	SU222	SU230	SU238	SU246	SU254	SU262	SU270	SU278	SU286
G	SU199	SU207	SU215	SU223	SU231	SU239	SU247	SU255	SU263	SU271	SU279	SU287
H	SU200	SU208	SU216	SU224	SU232	SU240	SU248	SU256	SU264	SU272	SU280	SU288

SN91096-4-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU289	SU297	SU305	SU313	SU321	SU329	SU337	SU345	SU353	SU361	SU369	SU377
B	SU290	SU298	SU306	SU314	SU322	SU330	SU338	SU346	SU354	SU362	SU370	SU378
C	SU291	SU299	SU307	SU315	SU323	SU331	SU339	SU347	SU355	SU363	SU371	SU379
D	SU292	SU300	SU308	SU316	SU324	SU332	SU340	SU348	SU356	SU364	SU372	SU380
E	SU293	SU301	SU309	SU317	SU325	SU333	SU341	SU349	SU357	SU365	SU373	SU381
F	SU294	SU302	SU310	SU318	SU326	SU334	SU342	SU350	SU358	SU366	SU374	SU382
G	SU295	SU303	SU311	SU319	SU327	SU335	SU343	SU351	SU359	SU367	SU375	SU383
H	SU296	SU304	SU312	SU320	SU328	SU336	SU344	SU352	SU360	SU368	SU376	SU384

SN91384-PLATES-B (All four plates; SN91096-5-PLATE, SN91096-6-PLATE, SN91096-7-PLATE, SN91096-8-PLATE are included in bundle).

SN91096-5-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU385	SU393	SU401	SU409	SU417	SU425	SU433	SU441	SU449	SU457	SU465	SU473
B	SU386	SU394	SU402	SU410	SU418	SU426	SU434	SU442	SU450	SU458	SU466	SU474
C	SU387	SU395	SU403	SU411	SU419	SU427	SU435	SU443	SU451	SU459	SU467	SU475
D	SU388	SU396	SU404	SU412	SU420	SU428	SU436	SU444	SU452	SU460	SU468	SU476
E	SU389	SU397	SU405	SU413	SU421	SU429	SU437	SU445	SU453	SU461	SU469	SU477
F	SU390	SU398	SU406	SU414	SU422	SU430	SU438	SU446	SU454	SU462	SU470	SU478
G	SU391	SU399	SU407	SU415	SU423	SU431	SU439	SU447	SU455	SU463	SU471	SU479
H	SU392	SU400	SU408	SU416	SU424	SU432	SU440	SU448	SU456	SU464	SU472	SU480

SN91096-6-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU481	SU489	SU497	SU505	SU513	SU521	SU529	SU537	SU545	SU553	SU561	SU569
B	SU482	SU490	SU498	SU506	SU514	SU522	SU530	SU538	SU546	SU554	SU562	SU570
C	SU483	SU491	SU499	SU507	SU515	SU523	SU531	SU539	SU547	SU555	SU563	SU571
D	SU484	SU492	SU500	SU508	SU516	SU524	SU532	SU540	SU548	SU556	SU564	SU572
E	SU485	SU493	SU501	SU509	SU517	SU525	SU533	SU541	SU549	SU557	SU565	SU573
F	SU486	SU494	SU502	SU510	SU518	SU526	SU534	SU542	SU550	SU558	SU566	SU574
G	SU487	SU495	SU503	SU511	SU519	SU527	SU535	SU543	SU551	SU559	SU567	SU575
H	SU488	SU496	SU504	SU512	SU520	SU528	SU536	SU544	SU552	SU560	SU568	SU576

SN91096-7-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU577	SU585	SU593	SU601	SU609	SU617	SU625	SU633	SU641	SU649	SU657	SU665
B	SU578	SU586	SU594	SU602	SU610	SU618	SU626	SU634	SU642	SU650	SU658	SU666
C	SU579	SU587	SU595	SU603	SU611	SU619	SU627	SU635	SU643	SU651	SU659	SU667
D	SU580	SU588	SU596	SU604	SU612	SU620	SU628	SU636	SU644	SU652	SU660	SU668
E	SU581	SU589	SU597	SU605	SU613	SU621	SU629	SU637	SU645	SU653	SU661	SU669
F	SU582	SU590	SU598	SU606	SU614	SU622	SU630	SU638	SU646	SU654	SU662	SU670
G	SU583	SU591	SU599	SU607	SU615	SU623	SU631	SU639	SU647	SU655	SU663	SU671
H	SU584	SU592	SU600	SU608	SU616	SU624	SU632	SU640	SU648	SU656	SU664	SU672

SN91096-8-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU673	SU681	SU689	SU697	SU705	SU713	SU721	SU729	SU737	SU745	SU753	SU761
B	SU674	SU682	SU690	SU698	SU706	SU714	SU722	SU730	SU738	SU746	SU754	SU762
C	SU675	SU683	SU691	SU699	SU707	SU715	SU723	SU731	SU739	SU747	SU755	SU763
D	SU676	SU684	SU692	SU700	SU708	SU716	SU724	SU732	SU740	SU748	SU756	SU764
E	SU677	SU685	SU693	SU701	SU709	SU717	SU725	SU733	SU741	SU749	SU757	SU765
F	SU678	SU686	SU694	SU702	SU710	SU718	SU726	SU734	SU742	SU750	SU758	SU766
G	SU679	SU687	SU695	SU703	SU711	SU719	SU727	SU735	SU743	SU751	SU759	SU767
H	SU680	SU688	SU696	SU704	SU712	SU720	SU728	SU736	SU744	SU752	SU760	SU768

SN91384-PLATES-C (All four plates; SN91096-9-PLATE, SN91096-10-PLATE, SN91096-11-PLATE, SN91096-12-PLATE are included in bundle).

SN91096-9-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU769	SU777	SU785	SU793	SU801	SU809	SU817	SU825	SU833	SU841	SU849	SU857
B	SU770	SU778	SU786	SU794	SU802	SU810	SU818	SU826	SU834	SU842	SU850	SU858
C	SU771	SU779	SU787	SU795	SU803	SU811	SU819	SU827	SU835	SU843	SU851	SU859
D	SU772	SU780	SU788	SU796	SU804	SU812	SU820	SU828	SU836	SU844	SU852	SU860
E	SU773	SU781	SU789	SU797	SU805	SU813	SU821	SU829	SU837	SU845	SU853	SU861
F	SU774	SU782	SU790	SU798	SU806	SU814	SU822	SU830	SU838	SU846	SU854	SU862
G	SU775	SU783	SU791	SU799	SU807	SU815	SU823	SU831	SU839	SU847	SU855	SU863
H	SU776	SU784	SU792	SU800	SU808	SU816	SU824	SU832	SU840	SU848	SU856	SU864

SN91096-10-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU865	SU873	SU881	SU889	SU897	SU905	SU913	SU921	SU929	SU937	SU945	SU953
B	SU866	SU874	SU882	SU890	SU898	SU906	SU914	SU922	SU930	SU938	SU946	SU954
C	SU867	SU875	SU883	SU891	SU899	SU907	SU915	SU923	SU931	SU939	SU947	SU955
D	SU868	SU876	SU884	SU892	SU900	SU908	SU916	SU924	SU932	SU940	SU948	SU956
E	SU869	SU877	SU885	SU893	SU901	SU909	SU917	SU925	SU933	SU941	SU949	SU957
F	SU870	SU878	SU886	SU894	SU902	SU910	SU918	SU926	SU934	SU942	SU950	SU958
G	SU871	SU879	SU887	SU895	SU903	SU911	SU919	SU927	SU935	SU943	SU951	SU959
H	SU872	SU880	SU888	SU896	SU904	SU912	SU920	SU928	SU936	SU944	SU952	SU960

SN91096-11-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU961	SU969	SU977	SU985	SU993	SU1001	SU1009	SU1017	SU1025	SU1033	SU1041	SU1049
B	SU962	SU970	SU978	SU986	SU994	SU1002	SU1010	SU1018	SU1026	SU1034	SU1042	SU1050
C	SU963	SU971	SU979	SU987	SU995	SU1003	SU1011	SU1019	SU1027	SU1035	SU1043	SU1051
D	SU964	SU972	SU980	SU988	SU996	SU1004	SU1012	SU1020	SU1028	SU1036	SU1044	SU1052
E	SU965	SU973	SU981	SU989	SU997	SU1005	SU1013	SU1021	SU1029	SU1037	SU1045	SU1053
F	SU966	SU974	SU982	SU990	SU998	SU1006	SU1014	SU1022	SU1030	SU1038	SU1046	SU1054
G	SU967	SU975	SU983	SU991	SU999	SU1007	SU1015	SU1023	SU1031	SU1039	SU1047	SU1055
H	SU968	SU976	SU984	SU992	SU1000	SU1008	SU1016	SU1024	SU1032	SU1040	SU1048	SU1056

SN91096-12-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU1057	SU1065	SU1073	SU1081	SU1089	SU1097	SU1105	SU1113	SU1121	SU1129	SU1137	SU1145
B	SU1058	SU1066	SU1074	SU1082	SU1090	SU1098	SU1106	SU1114	SU1122	SU1130	SU1138	SU1146
C	SU1059	SU1067	SU1075	SU1083	SU1091	SU1099	SU1107	SU1115	SU1123	SU1131	SU1139	SU1147
D	SU1060	SU1068	SU1076	SU1084	SU1092	SU1100	SU1108	SU1116	SU1124	SU1132	SU1140	SU1148
E	SU1061	SU1069	SU1077	SU1085	SU1093	SU1101	SU1109	SU1117	SU1125	SU1133	SU1141	SU1149
F	SU1062	SU1070	SU1078	SU1086	SU1094	SU1102	SU1110	SU1118	SU1126	SU1134	SU1142	SU1150
G	SU1063	SU1071	SU1079	SU1087	SU1095	SU1103	SU1111	SU1119	SU1127	SU1135	SU1143	SU1151
H	SU1064	SU1072	SU1080	SU1088	SU1096	SU1104	SU1112	SU1120	SU1128	SU1136	SU1144	SU1152

SN91384-PLATES-D (All four plates; SN91096-13-PLATE, SN91096-14-PLATE, SN91096-15-PLATE, SN91096-16-PLATE are included in bundle).

SN91096-13-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU1153	SU1161	SU1169	SU1177	SU1185	SU1193	SU1201	SU1209	SU1217	SU1225	SU1233	SU1241
B	SU1154	SU1162	SU1170	SU1178	SU1186	SU1194	SU1202	SU1210	SU1218	SU1226	SU1234	SU1242
C	SU1155	SU1163	SU1171	SU1179	SU1187	SU1195	SU1203	SU1211	SU1219	SU1227	SU1235	SU1243
D	SU1156	SU1164	SU1172	SU1180	SU1188	SU1196	SU1204	SU1212	SU1220	SU1228	SU1236	SU1244
E	SU1157	SU1165	SU1173	SU1181	SU1189	SU1197	SU1205	SU1213	SU1221	SU1229	SU1237	SU1245
F	SU1158	SU1166	SU1174	SU1182	SU1190	SU1198	SU1206	SU1214	SU1222	SU1230	SU1238	SU1246
G	SU1159	SU1167	SU1175	SU1183	SU1191	SU1199	SU1207	SU1215	SU1223	SU1231	SU1239	SU1247
H	SU1160	SU1168	SU1176	SU1184	SU1192	SU1200	SU1208	SU1216	SU1224	SU1232	SU1240	SU1248

SN91096-14-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU1249	SU1257	SU1265	SU1273	SU1281	SU1289	SU1297	SU1305	SU1313	SU1321	SU1329	SU1337
B	SU1250	SU1258	SU1266	SU1274	SU1282	SU1290	SU1298	SU1306	SU1314	SU1322	SU1330	SU1338
C	SU1251	SU1259	SU1267	SU1275	SU1283	SU1291	SU1299	SU1307	SU1315	SU1323	SU1331	SU1339
D	SU1252	SU1260	SU1268	SU1276	SU1284	SU1292	SU1300	SU1308	SU1316	SU1324	SU1332	SU1340
E	SU1253	SU1261	SU1269	SU1277	SU1285	SU1293	SU1301	SU1309	SU1317	SU1325	SU1333	SU1341
F	SU1254	SU1262	SU1270	SU1278	SU1286	SU1294	SU1302	SU1310	SU1318	SU1326	SU1334	SU1342
G	SU1255	SU1263	SU1271	SU1279	SU1287	SU1295	SU1303	SU1311	SU1319	SU1327	SU1335	SU1343
H	SU1256	SU1264	SU1272	SU1280	SU1288	SU1296	SU1304	SU1312	SU1320	SU1328	SU1336	SU1344

SN91096-15-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU1345	SU1353	SU1361	SU1369	SU1377	SU1385	SU1393	SU1401	SU1409	SU1417	SU1425	SU1433
B	SU1346	SU1354	SU1362	SU1370	SU1378	SU1386	SU1394	SU1402	SU1410	SU1418	SU1426	SU1434
C	SU1347	SU1355	SU1363	SU1371	SU1379	SU1387	SU1395	SU1403	SU1411	SU1419	SU1427	SU1435
D	SU1348	SU1356	SU1364	SU1372	SU1380	SU1388	SU1396	SU1404	SU1412	SU1420	SU1428	SU1436
E	SU1349	SU1357	SU1365	SU1373	SU1381	SU1389	SU1397	SU1405	SU1413	SU1421	SU1429	SU1437
F	SU1350	SU1358	SU1366	SU1374	SU1382	SU1390	SU1398	SU1406	SU1414	SU1422	SU1430	SU1438
G	SU1351	SU1359	SU1367	SU1375	SU1383	SU1391	SU1399	SU1407	SU1415	SU1423	SU1431	SU1439
H	SU1352	SU1360	SU1368	SU1376	SU1384	SU1392	SU1400	SU1408	SU1416	SU1424	SU1432	SU1440

SN91096-16-PLATE

	1	2	3	4	5	6	7	8	9	10	11	12
A	SU1441	SU1449	SU1457	SU1465	SU1473	SU1481	SU1489	SU1497	SU1505	SU1513	SU1521	SU1529
B	SU1442	SU1450	SU1458	SU1466	SU1474	SU1482	SU1490	SU1498	SU1506	SU1514	SU1522	SU1530
C	SU1443	SU1451	SU1459	SU1467	SU1475	SU1483	SU1491	SU1499	SU1507	SU1515	SU1523	SU1531
D	SU1444	SU1452	SU1460	SU1468	SU1476	SU1484	SU1492	SU1500	SU1508	SU1516	SU1524	SU1532
E	SU1445	SU1453	SU1461	SU1469	SU1477	SU1485	SU1493	SU1501	SU1509	SU1517	SU1525	SU1533
F	SU1446	SU1454	SU1462	SU1470	SU1478	SU1486	SU1494	SU1502	SU1510	SU1518	SU1526	SU1534
G	SU1447	SU1455	SU1463	SU1471	SU1479	SU1487	SU1495	SU1503	SU1511	SU1519	SU1527	SU1535
H	SU1448	SU1456	SU1464	SU1472	SU1480	SU1488	SU1496	SU1504	SU1512	SU1520	SU1528	SU1536

Section H: Product Ordering Information

Workflow Component	Product Name	Catalog Number
SNAP Core Kits	Swift Normalase Amplicon Panels (SNAP) Core Kit (96 rxns, no indexing)	SN-5X296
	Swift Normalase Amplicon Panels (SNAP) Core Kit (4x96 rxns Bundle, no indexing)	SN-5X384
Panel Specific Primers	SARS-CoV-2 Panel (primers only, 96 rxns)	COVG1-96
	SARS-CoV-2 Additional Genome Coverage Panel (primers only, 96 rxns)	COVG1V2-96
	SARS-CoV-2 S Gene Panel (primers only, 96 rxns)	SGENE-96
SNAP CDI Primers	SNAP Set 1A Combinatorial Dual Indexing Primers (96-plex, 96 rxns)	SN-5S1A96
	SNAP Set 1B Combinatorial Dual Indexing Primers (96-plex, 96 rxns)	SN-5S1B96
	SNAP Set 2A Combinatorial Dual Indexing Primers (96-plex, 96 rxns)	SN-5S2A96
	SNAP Set 2B Combinatorial Dual Indexing Primers (96-plex, 96 rxns)	SN-5S2B96
	SNAP Set S1AB-S2AB Combinatorial Dual Indexing Primers (384-plex, 4x96 rxns Bundle)	SN-5S0384
SNAP UDI Primer Plates	SNAP Unique Dual Indexing Primer Plate (96-plex, 96 rxns, SU001-SU096)	SN91096-1-PLATE
	SNAP Unique Dual Indexing Primer Plate (96-plex, 96 rxns, SU097-SU192)	SN91096-2-PLATE
	SNAP Unique Dual Indexing Primer Plate (96-plex, 96 rxns, SU193-SU288)	SN91096-3-PLATE
	SNAP Unique Dual Indexing Primer Plate (96-plex, 96 rxns, SU289-SU384)	SN91096-4-PLATE
	SNAP Unique Dual Indexing Primer Plates (384-plex, 4x96 rxns Bundle, SU001-SU384)	SN91384-PLATES
	SNAP Unique Dual Indexing Primer Plates (384-plex, 4x96 rxns Bundle, SU385-SU768)	SN91384-B-PLATES
	SNAP Unique Dual Indexing Primer Plates (384-plex, 4x96 rxns Bundle, SU769-SU1152)	SN91384-C-PLATES
	SNAP Unique Dual Indexing Primer Plates (384-plex, 4x96 rxns Bundle, SU1153-SU1536)	SN91384-D-PLATES

Support

For additional support, please contact Swift at TechSupport@swiftbio.com, or by phone: 734.330.2568 (9:00 am – 5:00 pm ET, Monday through Friday).

Revision History

Document #	Revision	Date	Description of Change
PRT-028	Version 1	5/20/2020	Initial release.
PRT-028	Version 2	7/10/2020	Amended Kit Contents, Introduction to Normalase, Calculator and Indexing tables.
PRT-028	Version 3	8/24/2020	Added panel COVG1V2-96 Cat No. and New Product Information.
PRT-028	Version 4	8/31/2020	Amended multiplex PCR annealing temperature and cleanup step 2 bead ratio; Added PCR cycle number option for low viral loads.
PRT-028	Version 5	10/28/2020	Addition of Swift 384 UDIs and new index information strategy
PRT-028	Version 6	01/18/2021	Addition of SARS-CoV-2 S Gene Panel and reformatting.
PRT-028	Version 7	02/10/2021	Corrected minor spelling errors.
PRT-028	Version 8	04/14/2021	Updated recommendations to sequencing depth.
PRT-028	Version 9	05/03/2021	Expanded indexes offering to 1536 UDIs. Updates to indexing layout plates and ordering information.

Notice to Purchaser: Limited License

For research use only. Not for use in diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations.

674 S. Wagner Road, Suite 100 • Ann Arbor, MI 48103 • 734.330.2568 • www.swiftbiosci.com

© 2021, Swift Biosciences, Inc. The Swift logo and Swift Normalase are trademarks of Swift Biosciences and are registered in the U.S. SNAP is covered by one or more claims of US Patent No. 10,316,359. This product is for Research Use Only. Illumina®, MiniSeq®, MiSeq®, NextSeq®, NovaSeq®, HiSeq®, and TruSeq® are registered trademarks of Illumina, Inc. SPRI®, SPRIPlate®, and SPRIselect® are trademarks of Beckman Coulter, Inc. NanoDrop®, Superscript®, and Qubit® and DynaMag® are registered trademarks of Thermo Fisher Scientific, Inc. Oligonucleotide adapter sequences are copyrighted © 2021 Illumina, Inc. All rights reserved. PRT-028 Rev 9.